


Dometic GmbH, In der Steinwiese 16, D-57074 Siegen

### **DS 300H, DS 600H Medical Refrigerator Hungary** **Minus degrees within the Medical Refrigerator**

If minus degrees at the above mentioned fridges occur, please check as follows:

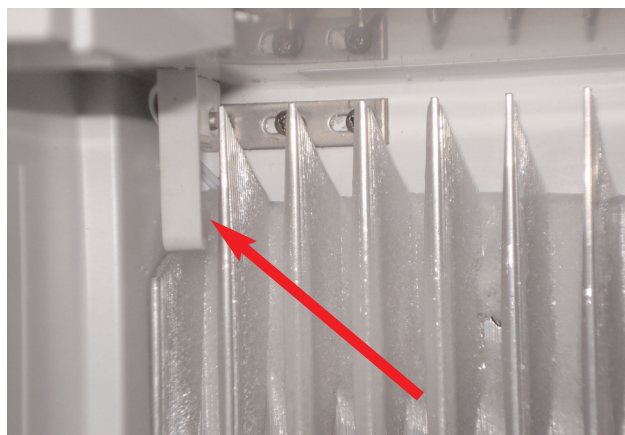

Icing at post evaporator.

Defrost the appliance in regular intervals.

The layer of frost should not come in contact with the temperature sensor. (Minimum distance 3mm)



The temperature sensor in the cooling compartment shall not be covered by the goods. The goods shall not be positioned straight at the backplane of the cooling compartment. Let air ventilate between the goods. (Minimum distance 5mm)

Please note, that a constant ventilation must be guaranteed if the fridge is built in. (Minimum free cross section 200cm<sup>2</sup>)